



PHYSICAL TESTER SOFTWARE


## EXPORT MANAGER

## Introducing simple and fast data export with the Erweka Export Manager.

The Erweka Export Manager allows users to seamlessly export test reports from devices such as SVM II, TAR II, EasyCheck, MultiCheck 6 and ZT 720/30 over their network and easily convert them to PDF and XML formats. The Windows software is installed on any PC within your laboratory network and automatically converts your test report data into PDF and XML files, doing away with the time-consuming manual scanning of printed reports.


But it doesn't stop there. Our user-friendly interface allows to define the desired destination folder, whether it's a local drive or a specific network location. By integration into the lab's own IT network, the security measures and data compliance protocols that apply to the network also apply to these folders, so that valuable data always remains protected. No compromise in data security or integrity: each exported file is given a specific name that includes the device type, serial number, date and time of data receipt.

Erweka Export Manager is a lightweight solution suitable for both new and existing devices. With its remarkable ability to receive data from multiple devices simultaneously, users save time, effort and money and streamline their laboratory workflows.

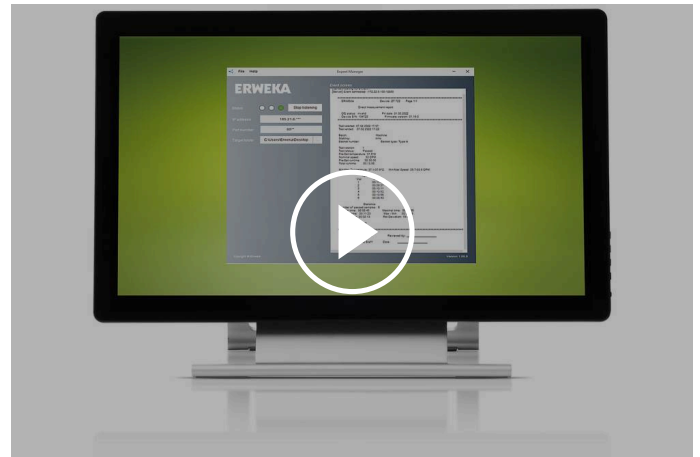

[Click to view video](#)

## EFFORTLESS SAVING ON PC OR IN NETWORK FOLDER STREAMLINING DATA COLLECTION FROM YOUR EQUIPMENT

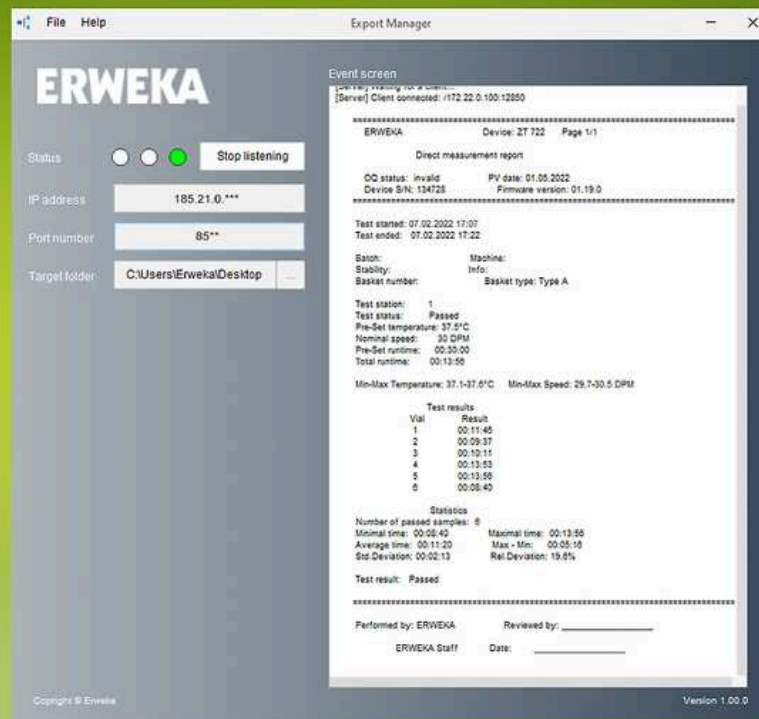

### STATUS

Shows the status of the network connection. Red: Values for "Port Number" and "Target Folder" are missing or incorrect. Yellow: The values entered are correct. Green: Data reception has been started. The software is ready to receive data from the device.

### START LISTENING

Switches the Export Manager to the "Waiting" status, in which data can be received. The button is active only after you have entered correct values for "Port Number" and "Target Folder".

## IP ADDRESS

Shows the IP address of the PC on which the Export Manager is running. It is determined automatically.

## PORT NUMBER

Sets the number of the port where data is received from the device.

## TARGET FOLDER

Sets the folder path where the protocol will be saved.  
Click to set a path.

# PHYSICAL TESTER SOFTWARE EXPORT MANAGER



The Export Manager software supports the following Erweka devices:

- | ZT 72x (from firmware version 1.20)
- | ZT 73x (from firmware version 2.02)
- | EasyCheck (from firmware version 1.38)
- | MultiCheck 6 (from firmware version 1.38)
- | SVM II (from firmware version 1.15)
- | TAR II (from firmware version 1.01)